



Doppler velocimetry

SPECIFICATION SHEET

AVAILABLE IN PXI

AVAILABLE IN MATRIQ

quantifiphotonics.com

FEATURES

Velocity measurement of fast-moving materials is essential to a growing number of scientific and technical fields including shock physics, the study of material properties, hypervelocity impact testing, detonics and shock wave experiments.

The Quantifi Photonics Doppler combines the key optical components for Photonic Doppler Velocimetry (PDV) measurements in a compact instrument to enable streamlined high channel count PDV test setups.



TARGET APPLICATIONS

- Study of material properties and impact response using shock physics.
- Replacement for Velocity Interferometer System for Any Reflector (VISAR) with an integrated fiber-based system.
- Measure velocities in dynamic experiments with high temporal precision.

A COMPLETE PDV MEASUREMENT SOLUTION

Combine the Doppler with the Laser and O2E to build a streamlined PDV test system. All you need to add is an optical probe and digitizer or oscilloscope with suitable bandwidth for the application.













LASER Tunable Laser Source

Available in PXI Available in MATRIQ

- Continuous Wave tunable laser source offering high-power output, narrow 100kHz linewidth and 0.01 pm resolution tunability.
- Whisper Mode is an optional feature to remotely disable frequency dither on each laser.

DOPPLER Photonic Doppler Velocimeter

Available in PXI Available in MATRIQ

- Available in standard and powercontrolled models.
- Power meter, variable optical attenuator and coupler all in one compact, space saving unit.

O2E Optical-to-Electrical Converter

Available in PXI Available in MATRIQ

- High-bandwidth, broadband optical-to-electrical converter with 1 or 2 channels, AC or DC coupling, and various conversion gain and operating wavelength ranges.
- Built-in amplifier eliminates the need for additional RF amplifier for significant cost savings per channel.

FIGURE 1

Figure 1 represents the configuration of the Doppler 1001 in PXI or MATRIQ form. It incorporates many of the key fiber optical components required to accurately condition the target and the reference laser and to mix the probe with the reference.


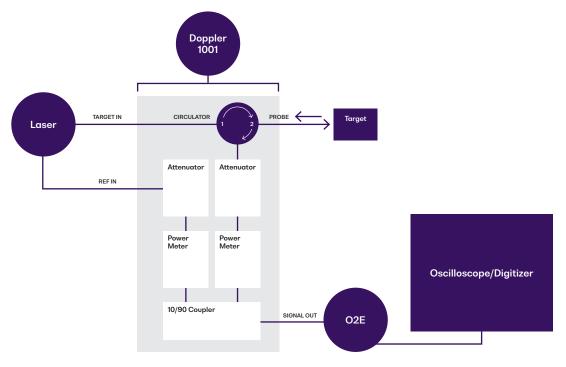


FIGURE 2

Figure 2 represents an alternative configuration of the Doppler instrument. The DOPPLER 1002 places the target input attenuator and power meter before the circulator. With the power control before the probe, this reduces the optical power transmitted to the probe and the DUT which may be desirable in some circumstances.

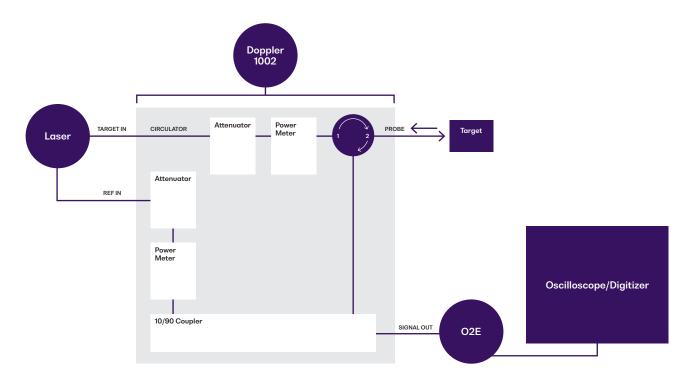

PXIe - MODULAR

Our expanding range of PXIe optical test solutions are used by customers in mixed-signal test and measurement systems, reducing complexity, lowering the cost of test and accelerating time to market.

- Multi vendor, open standard with over 2500 PXI modules available
- Advanced timing and synchronization capabilities across instruments
- Low latency, high performance processing and fast data throughput
- Design and build scalable, high channel count systems
- Small footprint and lower power consumption



MATRIQ - COMPACT & PORTABLE

The MATRIQ series provides the same high-performance test capabilities of our PXIe modules in an compact benchtop design. MATRIQ instruments are simple to setup and easy to operate, making them the perfect choice for your optical lab or test bench.


- Same performance and control as our PXIe modules
- Plug and play with USB or Ethernet connectivity
- Control via the web-based GUI, COHESIONUI or SCPI commands
- Compact and portable design saves benchtop space




DOPPLER TECHNICAL SPECIFICATIONS



MATRIQ - COMPACT & PORTABLE

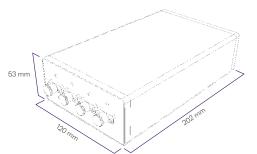

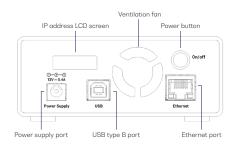





DOPPLER-1001-1-FC-MTRQ

PXI - 8-CHANNEL PDV SETUP

The below shows an example configuration of an 8-channel Doppler setup. It consists of four Laser 1002 laser sources (Target and Reference), and eight Doppler 1001 modules. The output of the Doppler 1001 module is sent to a four high channel and high bandwidth O2E 1101 where the probe and reference laser beat. The electrical beat signal is then digitized with an oscilloscope.

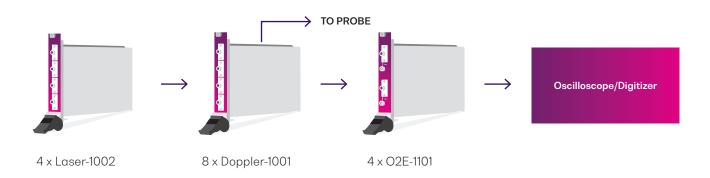

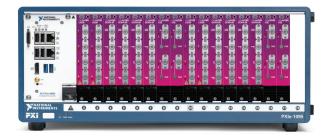

8-channel PDV system in a single 18-slot chassis

DOPPLER TECHNICAL SPECIFICATIONS

General Specifications	PXI	MATRIQ	
Bus connection	PXIe	USB and Ethernet	
Number of PXI slots	1	-	
Fiber type	SMF-28	SMF-28	
Optical connector type	FC/PC,SC/PC,FC/APC or SC/APC	FC/PC,SC/PC,FC/APC or SC/APC	
Dimensions H x W x D	130 x 20 x 215 mm 5.1 x 0.8 x 8.5 inches	53 x 120 x 202 mm 2.1 x 4.7 x 8.0 inches	
Weight	~ 1 kg ~ 2.2 lbs	~ 1.1 kg ~ 2.4 lbs	
Operating temperature range	5 °C to 45 °C 41 °F to 113 °F		
Storage temperature range	-40 °C to 70 °C -40 °F to 158 °F		

Model Number	1001	1002	1001	1002
Wavelength	1520 to 1650 nm		1520 to 1650 nm	
Probe output operating range	-50 to +20 dBm		-50 to +20 dBm	
Reference output operating range	-60 to +10 dBm		-60 to +10 dBm	
Damage level	+23 dBm		+23 dBm	
Insertion loss ³ target input to probe output	< 3 dB	< 5 dB	< 3 dB	< 5 dB
Insertion loss ³ probe input to output	< 5 dB	< 3 dB	< 5 dB	< 3 dB
Insertion loss ³ reference input to output	< 14 dB	< 14 dB	< 14 dB	< 14 dB
Wavelength dependent loss	< 0.02 dB/nm		< 0.02 dB/nm	
Return loss ³	> 45 dB		> 45 dB	
Circulator Directivity	> 45 dB		> 45 dB	
Warm-up time	< 20 minutes		< 20 minutes	

Attenuator	1001	1002	1001	1002
Calibration wavelength	1550 nm		1550 nm	
Attenuation range typical ⁵	> 46 dB		> 46 dB	
Attenuation range guaranteed ⁵	> 40 dB		> 40 dB	
Resolution	0.01 dB		0.01 dB	
Attenuation speed	0.1 to 1000 dB/s		0.1 to 1000 dB/s	


Power Meter	1001	1002	1001	1002
Calibration wavelength	1550 nm		1550 nm	
Polarization dependent responsivity ^{2,3}	< 0.2 dB		< 0.2 dB	
Linearity ^{2,5}	± 0.1 dB		<u>+</u> 0.1 dB	
Total uncertainty ^{23,5}	<u>+</u> 0.34 dB (Typical); <u>+</u> 0.55 dB (Max)		<u>+</u> 0.34 dB (Typical); <u>+</u> 0.55 dB (Max)	
Averaging time	100 µs to 10 s		100 µs to 10 s	
Resolution	0.01 dB		0.01 dB	
Data logging	1 to 1024 per channel		1 to 1024 per channel	

 Notes

 1.
 Specifications are valid at 23 °C ± 3 °C

 2.
 + 10dBm to -40dBm, 23 °C

Excluding connectors
 At calibration wavelengths

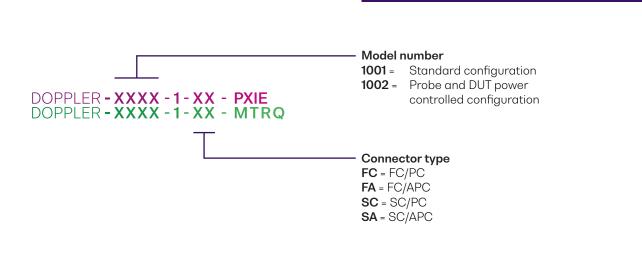

WARRANTY INFORMATION

ORDERING INFORMATION

This product comes with a standard 1 year warranty.

With an **extended warranty and calibration plan** you'll spend more time focused on your priorities and less time worrying about maintenance.

Your choice: add a **3 or 5 year extended** warranty when you buy.



Guarantee performance

Ensure your equipment is operating at the best it can be for reliable and accurate results.

Lower cost of ownership

Lock in savings and maximise your testing budget with a lower base cost of ownership.

Peace of mind

Spend less time worrying about maintenance and more on generating results.

CALIBRATION PLANS FOR ADDITIONAL DISCOUNTS

Order a **calibration plan** when purchasing your Quantifi Photonics instruments and get additional discounts.

10% Discount

On calibrations ordered at the time of purchase.

25% Discount

Add on an extended warranty and receive a 25% discount on calibrations.

Over time and with regular use, all optical parts and connectors require re-calibration and maintenance to guarantee accurate and reliable performance. We recommend Quantifi Photonics optical instruments are re-calibrated every 12 months. With an instrument calibration performed by Quantifi Photonics technicians you receive:

- Comprehensive calibration to factory specifications
- End-to-end inspection to ensure all instrument functions are working and connectors are clean
- Firmware, software and documentation updates
- Certificate of calibration which includes detailed test results

How to do I secure my extended warranty or calibration plan?

Contact your Quantifi Photonics sales representative or email sales@quantifiphotonics.com

Extended warranties and calibration plans must be ordered at the time of purchase and are available only for Quantifi Photonics' products. The 25% calibration discount only applies to calibrations while the product is covered by the extended warranty period.

CATALOGUE

Our portfolio of optical & electro-optical test modules is rapidly expanding to meet a wide range of customer requirements and applications.

Tunable Laser Sources

Versatile telecom laser sources with full tunability across C or L bands. Narrow 100 kHz linewidth, up to 16.5 dBm of power, optional whisper mode to disable frequency dither.

Superluminescent Diode **Broadband Light Source**

Super-luminescent LED light source with high output power, large bandwidth and low spectral ripple and various wavelenaths.

Polarization Controller & Scrambler

High-speed automated polarization control with broad . wavelength coverage from 1260nm to 1650nm, low insertion loss and back reflection. Full remote control via intuitive GUI LabVIEW or SCPI.

Digital Sampling Oscilloscope (DSO)

Digital equivalent-time sampling oscilloscope (DSO) with high-quality precision timebase and low jitter mode, available in 1 or 2 channels in a compact benchtop instrument.

Photonic Doppler Velocimeter (PDV)

Purpose-built module for Photonic Doppler Velocimetry (PDV). A circulator, two VOAs and a passive coupler all built into one compact module.

Fixed Wavelength Laser Sources

Highly customizable laser platform. Select required wavelength, power and fiber type for a customized solution.

Optical-to-Electrical Converter

High bandwidth, broadband O-to-E converter. Available in a range of configurations; choose from 1 or 2 channels, AC or DC coupling and various conversion agin and operating wavelength ranges.

Optical Power Meters

Fast terminating or inline monitoring of optical signal power from -60 to +10 dBm across 750 – 1700 nm wavelengths. Model with logarithmic analog output for applications such as silicon photonics fiber alignment.

Bit Error Rate Tester (BERT)

4 or 8-channel Pulse Pattern Generator and Error Detector at rates up to 29 Gbps for the design, characterization and production of optical transceivers and optoelectrical components.

Optical Switch

Proven reliability and fast switching time. Wide variety of switch onfigurations: 1x4, 1x16, 16x16 and more. Models support SMF, MMF and PMF.



Swept, Tunable Continuous Wave Laser

Swept, tunable continuous wave (CW) laser source with 0.01 dB power stability and 400 nm/s high-speed scan rate for R&D and production testing





Fast attenuation speed with low insertion loss and built-in power monitoring. Operates in fixed attenuation or constant output power modes. Models support SMF, MMF and PMF connector types.



Cost-effective, spectral measurement in a compact module with built-in analysis for: SMSR, OSNR & spectral width. Targeted wavelengths for specific applications in O band, C band & L band.

Passive Component Integration

Integrate passive optical components of your choice such as WDM couplers, splitters, band-pass filters, PM beamsplitters and circulators. Models support SMF, MMF and PMF.

Passive Component Storage

Protect and store your own passive fiber optic components such as splitters, connector adaptor patchcords, WDM couplers, and isolators in one







handy module.



For more details visit quantifiphotonics.com/products















Test. Measure. Solve[™]

Quantifi Photonics is transforming the world of photonics test and measurement. Our portfolio of optical and electrical test instruments is rapidly expanding to meet the needs of engineers and scientists around the globe. From enabling ground-breaking experiments to driving highly efficient production testing, you'll find us working with customers to solve complex problems with experience and innovation.

To find out more, get in touch with us today.

General Enquiries Technical Support Phone - NZ Phone - USA sales@quantifiphotonics.com support@quantifiphotonics.com +64 9 478 4849 +1-800-803-8872



quantifiphotonics.com

Quantifi Photonics Ltd © 2024. All rights reserved. No part of this publication may be reproduced, adapted, or translated in any form or by any means without the prior permission from Quantifi Photonics Ltd. All specifications are subject to change without notice. Please contact Quantifi Photonics for the latest information.